

## **Optojump Power-Up Guide**

- Turn on power bar, laptop, and the Optojump bars
- Plug in USB/Serial adaptor
- On laptop, click Start
- Click Settings if available then Control Panel. Or just open the Control Panel
- Click System
- Click Hardware tab
- Open Device Manager
- Expand Ports (COM+LPT); verify which COM port the USB/Serial adaptor is plugged into
- Close all of these windows
- Open Optojump software on the Desktop
- Click on Configuration tab
- Click Serial Port (when window opens click on box for correct COM port being used)
- Click OK
- Click on Configuration tab
- Click Optojump Test
- Enter 1 for Number of Bars
- Click Enter
- Press Start
- You must see this message: 'Trial successfully terminated' - (if there is an error, check for dust/debris in grid area and run test again)
- Click Exit
- Optojump is now ready for use